



International Filter Products

Revolutionary systems for your industry

Sterilization & Disinfection

SteriPure-G can maintain integrity after being exposed to a variety of harsh sterilization methods, for results see table below.

Autoclave	124 °C, 30 minutes, 100 Cycles
Steam in Place	124 °C, 30 minutes, 100 Cycles
Hot Water	85 °C, 60 minutes, 80 Cycles
NaOH Solution	85 °C, 3.0%, 80 Hours
HNO ₃	25 °C, 0.5, 100 Hours
H ₃ PO ₄	25 °C, 0.5, 100 Hours
H ₂ O ₂	25 °C, 1,200 ppm, 500 Hours
KHSO ₄ Solution	25 °C, 1,600 ppm, 500 Hours

The common sterilization and disinfection methods above have been tested under controlled laboratory conditions.

Note: This Cartridge has a glass fiber pre-filter laid over the PES membrane final filter and both layers are pleated together.

7914 Ajay Dr., Sun Valley, CA 91352
www.internationalfilterproducts.com
sales@internationalfilterproducts.com
Phone: (818) 841-2702
Fax: (818) 841-2947