STERILE CAPSULE FILTER

These filters are designed for applications where sterility of the affluent is mandatory (Sterile Compounding and Pharmaceutical use). The product contains a pharmaceutical grade PTFE membrane (Sterilizing Grade) and the filters are built in a 10,000-class clean room, flushed, dried and 100% integrity tested prior to packaging. The product is factory sterilized by Ethylene Oxide gas (ETO) ensuring sterility.

Technical Data Sheet

Catalog Number: SKLF100020H2H2H-PH-ETO & SKLF100020HLF2H-PH-ETO & SKLF100020HLFLM-PH-ETO







Materials of Construction

Filter Membrane: 1.0 um PTFE Pre-filter / 0.2 um final Pharmaceutical Grade PTFE Sterilizing Grade - Hydrophobic

Supports: Polypropylene

Core: Polypropylene

Capsule Body: Polypropylene

Filter Dimensions and Specifications

Outer Diameter: 71.5 mm (2.81")

Length, Fitting to Fitting: 115 mm (4.52") (2H2H)/ 102 mm (4.01") (LF2H)

Filter Area: 680 cm² (105.4 in²)

Sterilization: The filter is factory sterilized by Ethylene Oxide Gas (ETO)


Inlet Fitting: 2H2H: ¼" – ½" Tapered Hose Barb; **LF2H:** - Luer Loc Female

Outlet Fitting: ¼" – ½" Tapered Hose Barb

Volume Filtered: 15 L to 33.0 L





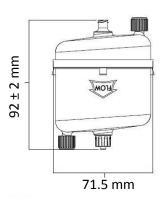

Integrity Data

| Pore Size | Min, Bubble Point, 25°C | Challenge Microorganisms | CFU per cm ² |
|-----------|-------------------------|--------------------------|-------------------------|
| 0.20 μm | ≥3.5 bar (50 psi) | Brevundimonas Diminuta | 7 |
| | | (ATCC 19146) | |

Operating Parameters

Maximum Working Pressure: 5.5 bar (80 Psi) @ 25 °C (77 °F)

Maximum differential Pressure (forward): 4.1 bar (60 Psi) @ 25 °C (77 °F)

Maximum differential Pressure (reverse): 2.1 bar (30 Psi) @ 25 °C (77 °F)

Maximum Operating Temperature: 80 °C (176 °F)

Recommended Replacement Pressure: 2.4 bar (35.3 Psi)

Safety

UPS Bacteria Endotoxins: ≤25 EU/ml

Factory Sterilized: Ethylene Oxide Gas (ETO)

Product was 100% Integrity Tested and flushed with filtered water @ 0.05~um

Regulatory

Complies with USP 797 Guidelines

ASTM F838-05 - Bacterial Retention

FDA 21 CFR 177.1655

USP Class VI - Biological Reactivity

ISO 10993-Part 1.5 - Cytotoxicity

ISO 14001; ISO 13485; OHSMS 18001 Certified.

ASTM Hemolysis testing



www.internationalfilterproducts.com 11070 Fleetwood St. Unit B Sun Valley, CA 91352 Tel: (818) 504-8115 - Office

Tel: (818) 917-9538 - Cell

E-mail: sales@internationalfilterproducts.com